

## Label printer settings

Brother QL-Serie



### Installing the drivers

First install the drivers. You will find these on the USB flash drive supplied along with the device under *Useful items/Brother drivers*. Select your model and start the application D\_SETUP.exe. Follow the installation instructions. Alternatively, you can find the latest drivers online at the [Brother Solution Centre](#).

### Setting up in the Apo-Ident software

If you have installed the drivers successfully, you can now choose your printer from the **Standard label printer** list (Brother QL-700 or older models) under **Label Printer Settings**.

### Continuous Paper Settings DK-22205

Choose the following settings:

- Page size: 62mm
- Orientation: Rotated by 0°

Advanced Layout Settings:

- Label width / mm: 62,0
- Label height / mm: 35,0
- X-offset / mm: 0,0
- Y-offset / mm: 0,0
- Scaling factor: 1,00





Now click on the left tool icon „Open printer settings“. Change the following settings in the dialogue window that opens:

- Paper Size: 62mm
- Length: 35.0
- Belt feed: 3.0
- Alignment: Portrait format
- Quality: Prioritise print quality 300 x 300 dpi

Click first on **<Apply>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

**Note:** You can check your settings by starting a test print. To do this, click on the middle icon "Print test label".

If your test print was successful, click **<Close>**. Your settings are accepted and saved.

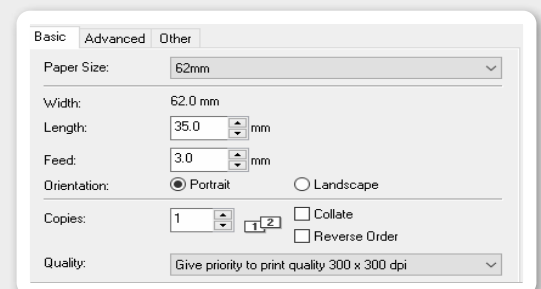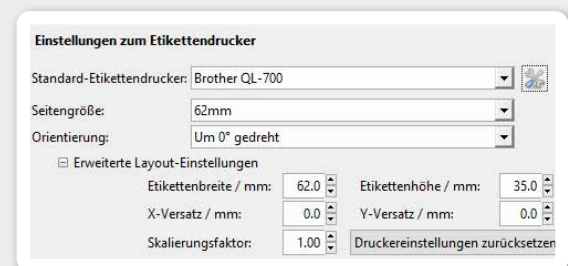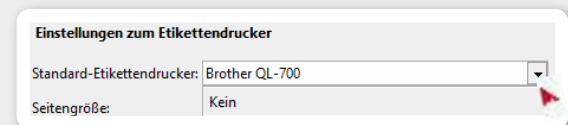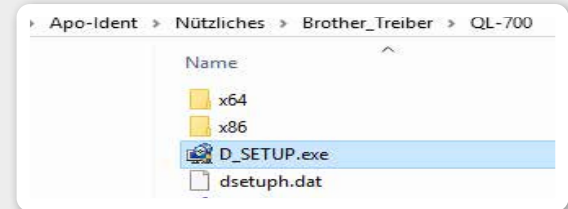

## Label printer settings

Brother QL-Serie



### Settings for single labels DK-11201

Choose the following settings:

- Page Size: 29 mm x 90 mm
- Orientation: Rotated by 90°

Advanced Layout Settings:

- Label width / mm: 29.0
- Label height / mm: 89.9
- X-offset / mm: 0.0
- Y-offset / mm: 0.0
- Scaling factor: 1.00



Now click on the left **tool icon** „Open printer settings“. In the dialogue window that opens, modify the following settings:

- Paper size: 29 mm x 90 mm
- Alignment: Portrait format
- Quality: Prioritise print quality 300 x 300 dpi

Click first on **<Apply>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

**Note:** You can check your settings by starting a test print. To do this, click on the middle icon „Print test label“.

If your test print was successful, click **<Close>**. Your settings are accepted and saved.

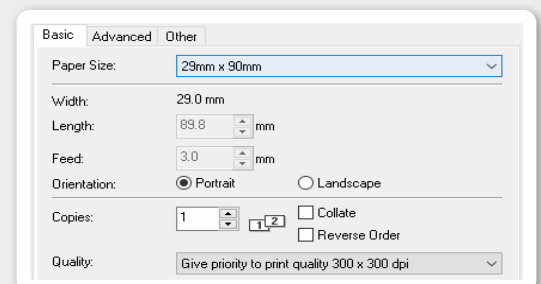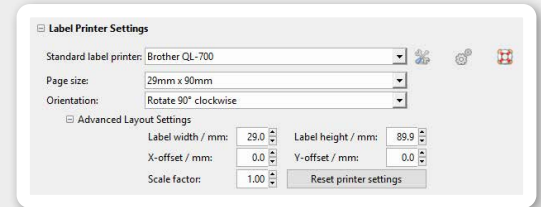

## Label printer settings

DYMO LabelWriter 450 for single labels



### Installing the driver

First install the driver. You will find these online at the [DYMO Support Center](#). After installing the printer driver, connect the printer to your PC.

### Setting up in the Apo-Ident software

If you have installed the drivers successfully, you can now choose your printer from the default label printer list under **<Label Printer Settings>**.

### Settings for single labels 99012

Choose the following settings:

- Default label printer: DYMO LabelWriter 450
- Page Size: 99012 Large Address
- Orientation: Rotate 0°

Advanced Layout Settings:

- Label width / mm: 35.8
- Label height / mm: 88.4
- X-offset / mm: 0.0
- Y-offset / mm: 0.0

Scaling factor: 2.20

Now click on the left tool icon **<Open printer settings>**. Change the following settings in the dialogue window that opens:

- Orientation: Landscape
- Page Order: Front to back




Click **<Advanced>** to make the following setting:

- Paper/Output: 99012 Large Address

Click first on **<OK>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

You can check your settings by starting a test print. To do this, click on the middle icon **<Print test label>**. If your test print was successful, click **<Close>**. Your settings are accepted and saved.

**Note:** These instructions only apply to the label printer DYMO LabelWriter 450 with labels 99012. With other DYMO models (e.g. Turbo, Twin Turbo, etc.) the label settings may differ.

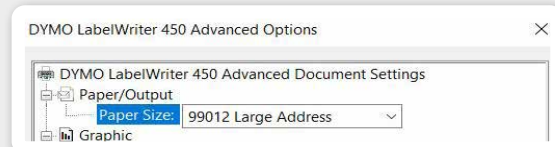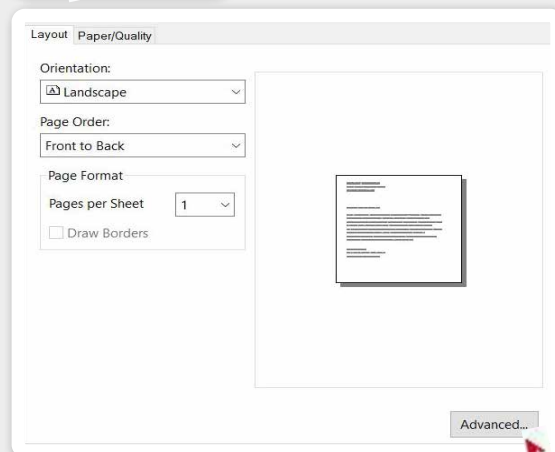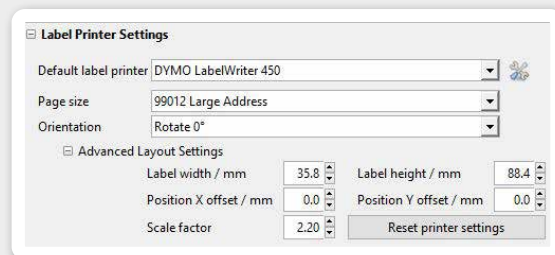

Apo-Ident ist eine Marke der  
HiperScan GmbH



HiperScan GmbH  
Weißeritzstraße 3  
01067 Dresden  
Germany

Phone: +49 351 212496-33  
Fax: +49 351 212496-99  
info@apo-ident.de  
www.apo-ident.de