

Label-printer settings

Installing driver software

First install the drivers, which can be found under Useful/Brother drivers on the USB stick supplied. Select your model and launch the D_SETUP.exe application. Follow the installation instructions. Alternatively, the latest drivers can also be found online in the [Brother Solution Centre](#).

Setting up the Apo-Ident software

Once you have successfully installed the drivers, you can select your printer from the list of **Standard label printers** (Brother QL-700 or older models) under **Label-printer settings**.

Settings with continuous DK-22205

Select the following settings:

- Page size: 62 mm
- Orientation: Rotate 0°

Advanced Layout settings

- Label width / mm: 62.0
- Label height / mm: 35.0
- X offset / mm: 0.0
- Y offset / mm: 0.0
- Scale factor: 1.00





Click the 'Open printer settings' **tool icon** on the left. Enter the following data in the dialogue box that appears:

- Paper size: 62 mm
- Length: 35.0
- Feed: 3.0
- Orientation: Portrait
- Quality: Give priority to print quality 300 x 300 dpi

First click **<Apply>** and confirm by clicking **<OK>**. This will take you back to the Apo-Ident software settings.

Note: You can check your settings by starting a test print. To do so, click the 'Print test label' icon in the middle.

If your test print was successful, click **<Close>**. Your settings will now be applied and saved.

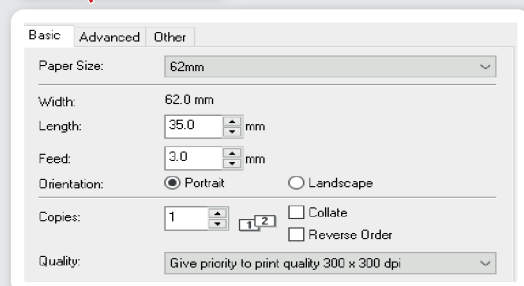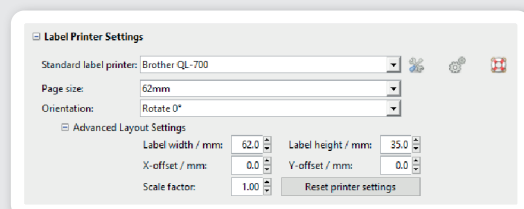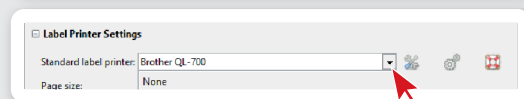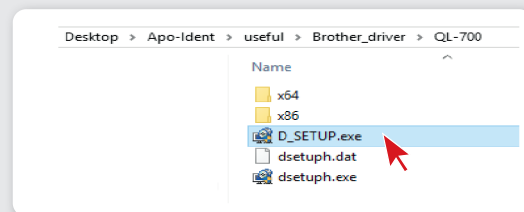

Settings with single labels DK-11201

Select the following settings:

- Page size: 29 mm x 90 mm
- Orientation: Rotate 90° clockwise

Advanced Layout settings

- Label width / mm: 29.0
- Label height / mm: 89.9
- X offset / mm: 0.0
- Y offset / mm: 0.0
- Scale factor: 1.00

Click the 'Open printer settings' **tool icon** on the left. Enter the following data in the dialogue box that appears:

- Paper size: 29 mm x 90 mm
- Orientation: Portrait
- Quality: Give priority to print quality 300 x 300 dpi

First click **<Apply>** and confirm by clicking **<OK>**. This will take you back to the Apo-Ident software settings.

Note: You can check your settings by starting a test print. To do so, click the 'Print test label' icon in the middle.

If your test print was successful, click **<Close>**. Your settings will now be applied and saved.

