

Label printer settings

Brother QL-Serie



Installing the drivers

First install the drivers. You will find these on the USB flash drive supplied along with the device under *Useful items/Brother drivers*. Select your model and start the application D_SETUP.exe. Follow the installation instructions. Alternatively, you can find the latest drivers online at the [Brother Solution Centre](#).

Setting up in the Apo-Ident software

If you have installed the drivers successfully, you can now choose your printer from the **Standard label printer** list (Brother QL-700 or older models) under **Label Printer Settings**.

Continuous Paper Settings DK-22205

Choose the following settings:

- Page size: 62mm
- Orientation: Rotated by 0°

Advanced Layout Settings:

- Label width / mm: 62,0
- Label height / mm: 35,0
- X-offset / mm: 0,0
- Y-offset / mm: 0,0
- Scaling factor: 1,00





Now click on the left tool icon „Open printer settings“. Change the following settings in the dialogue window that opens:

- Paper Size: 62mm
- Length: 35.0
- Belt feed: 3.0
- Alignment: Portrait format
- Quality: Prioritise print quality 300 x 300 dpi

Click first on **<Apply>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

Note: You can check your settings by starting a test print. To do this, click on the middle icon "Print test label".

If your test print was successful, click **<Close>**. Your settings are accepted and saved.

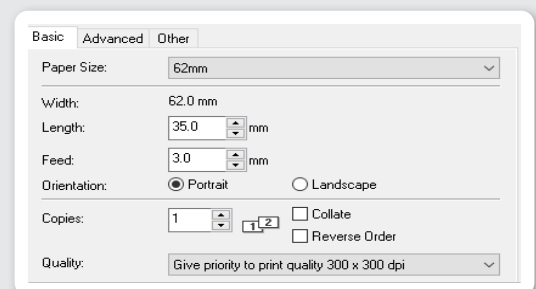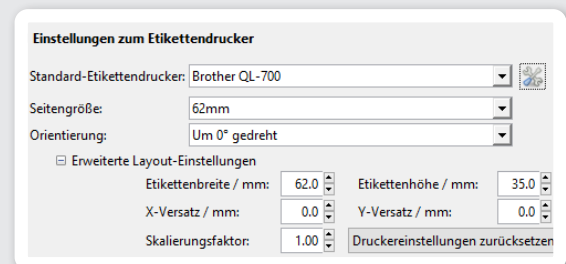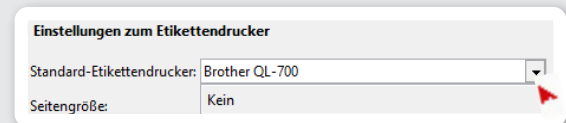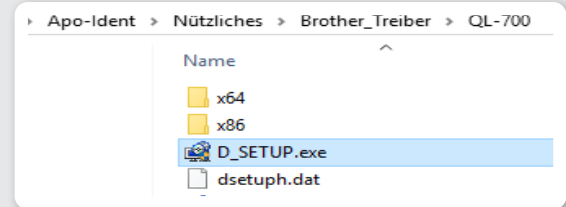

Label printer settings

Brother QL-Serie



Settings for single labels DK-11201

Choose the following settings:

- Page Size: 29 mm x 90 mm
- Orientation: Rotated by 90°

Advanced Layout Settings:

- Label width / mm: 29.0
- Label height / mm: 89.9
- X-offset / mm: 0.0
- Y-offset / mm: 0.0
- Scaling factor: 1.00



Now click on the left **tool icon** „Open printer settings“. In the dialogue window that opens, modify the following settings:

- Paper size: 29 mm x 90 mm
- Alignment: Portrait format
- Quality: Prioritise print quality 300 x 300 dpi

Click first on **<Apply>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

Note: You can check your settings by starting a test print. To do this, click on the middle icon „Print test label“.

If your test print was successful, click **<Close>**. Your settings are accepted and saved.

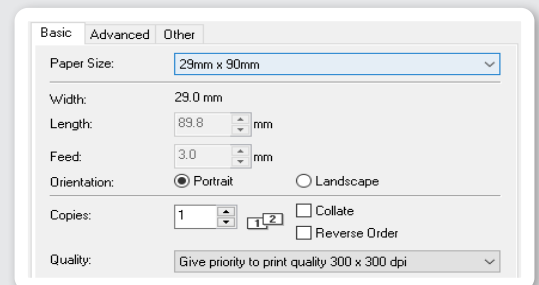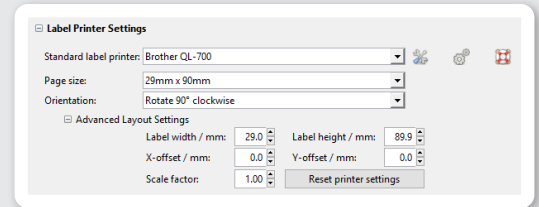

Label printer settings

DYMO LabelWriter 450 and 550 for single labels



Installing the driver

First install the driver. You will find these online at the [DYMO Support Center](#). After installing the printer driver, connect the printer to your PC.

Setting up in the Apo-Ident software

If you have installed the drivers successfully, you can now choose your printer from the default label printer list under **<Label Printer Settings>**.

Settings for single labels 99012

Choose the following settings:

- Default label printer: DYMO LabelWriter 450 **or** DYMO LabelWriter 550
- Page Size: 99012 Large Address
- Orientation: Rotate 0°

Advanced Layout Settings:

- Label width / mm: 35.8
- Label height / mm: 88.4
- X-offset / mm: 0.0
- Y-offset / mm: 0.0

Scaling factor: 2.20

Now click on the left tool icon **<Open printer settings>**. Change the following settings in the dialogue window that opens:

- Orientation: Landscape
- Page Order: Front to back




Click **<Advanced>** to make the following setting:

- Paper/Output: 99012 Large Address

Click first on **<OK>** and then confirm with **<OK>**. You are now back in the settings of the Apo-Ident software.

You can check your settings by starting a test print. To do this, click on the middle icon **<Print test label>**. If your test print was successful, click **<Close>**. Your settings are accepted and saved.

Note: These instructions only apply to the label printer DYMO LabelWriter 450/550 with labels 99012. With other DYMO models (e.g. Turbo, Twin Turbo, etc.) the label settings may differ.

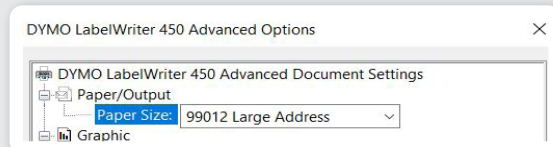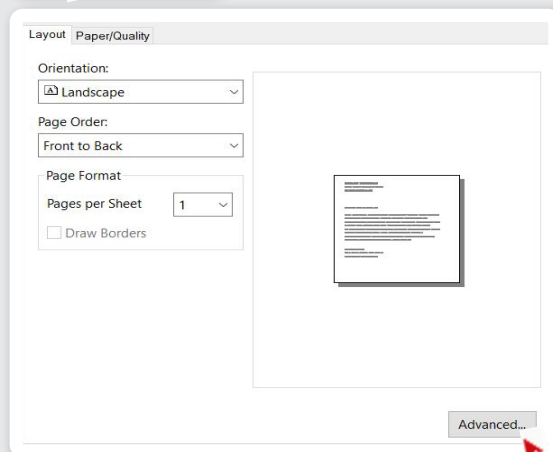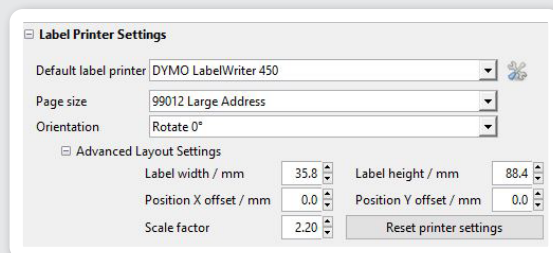

Apo-Ident is a trademark of
HiperScan GmbH



HiperScan GmbH
Gerokstraße 13
01307 Dresden
Germany

Phone: +49 351 212496-33
Fax: +49 351 212496-99
info@apo-ident.de
www.apo-ident.de